

CellShip®

A fully defined, xeno-free cell transportation medium
A simple and low-cost way to transport or store cells at ambient temperatures - your viable alternative to cryopreservation.


LSP
LIFE SCIENCE PRODUCTION

FEATURES

- A 'ready-to-use' transport and short-term storage medium, suitable for a wide range of cells
- Provides flexibility to highly time critical processes
- Removes the cost and logistic challenges associated with shipment on dry ice
- Provides stable pH-buffering of pH 7.20- 7.45 for temperatures between +2°C to +37°C
- Xeno-free and DMSO-free

Distributed by:

CliniSciences Group

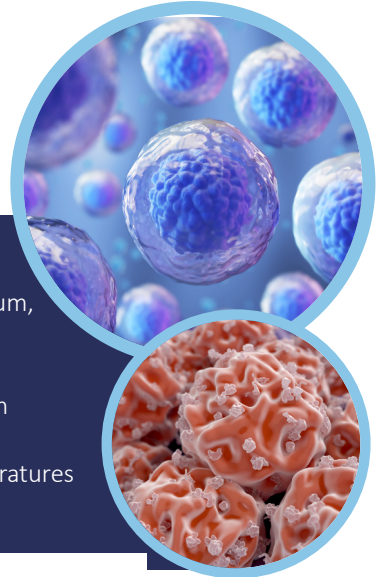

The protocol requires minimal cell manipulation, reducing the risk of contamination and improving efficiency within the laboratory.

Transport

Culture cells using standard growth media and conditions.

Harvest cells using your standard method for sub-culturing and centrifuge to remove supernatant.

Wash cells once in CellShip®.

Resuspend cell pellets in 2 mL of CellShip® and transfer to a 2 mL cryotube for transport.

Recovery

Gently resuspended the cells and transfer to the appropriate volume of complete medium and culture using standard conditions.

CellShip does not need to be removed before recovery!



Ordering Information

SCS-002X	CellShip Cell Transport Medium with Insulin 25mL, 50mL and 100mL
SCS-004X	CellShip Cell Transport Medium without Insulin 25mL, 50mL and 100mL

THIS PRODUCT IS CURRENTLY SOLD FOR RESEARCH PURPOSES ONLY

NOTE: the 'X' in the product code denotes pack size. Please state pack size required at time of order. Please note there is a charge per delivery.

Distributed by:

CliniSciences Group

CliniSciences Group

Netherlands

Company: CliniSciences B.V.
Address: Krijnburgstraat 137A,
1016RG Amsterdam, Netherlands
Telephone: +31 85 2082 351
Fax: +31 85 2082 353
Email: netherlands@clinisciences.com
Web: <http://www.clinisciences.com>



Norway

Company: CliniSciences ApS
Address: Østertorvegade 220, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 084
Email: norpo@clinisciences.com
Web: <http://www.clinisciences.com>



Poland

Company: CliniSciences sp.Z.o.o.
Address: ul. Roślinnicza Włódka Piłcockiego 67
lok. 200 - 02-701 Warszawa - Poland
Telephone: +48 22 307 0535
Fax: +48 22 307 0532
Email: polaka@clinisciences.com
Web: <http://www.clinisciences.com>



Austria

Company: CliniSciences GmbH
Address: Stammersgasse 76, A-1160
Wien - Austria
Telephone: +43 720 115 680
Fax: +43 720 115 577
Email: austorro@clinisciences.com
Web: <http://www.clinisciences.com>



Belgium

Company: CliniSciences S.R.L.
Address: Avenue Stalingrad 52, 1000
Brussels - Belgium
Telephone: +32 2 31 60 800
Fax: +32 2 31 60 801
Email: belgium@clinisciences.com
Web: <http://www.clinisciences.com>



Portugal

Company: Quimigén Unipessoal LDA
Address: Rua Almeida Negreiros, Lote 6, Loja 14,
2015-275 Alverca Do Ribatejo - Portugal
Telephone: +351 30 8808 050
Fax: +351 30 8808 052
Email: info@quimigen.com
Web: <http://www.quimigen.pt>



Spain

Company: CliniSciences Lab Soluções
Address: C/ Herranias del Moral 13
(Baño E), 28019, Madrid - Spain
Telephone: +34 91 269 40 65
Fax: +34 91 269 40 74
Email: spanna@clinisciences.com
Web: <http://www.clinisciences.com>



Sweden

Company: CliniSciences ApS
Address: Østertorvegade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: swedpo@clinisciences.com
Web: <http://www.clinisciences.com>



Finland

Company: CliniSciences ApS
Address: Østertorvegade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: suomi@clinisciences.com
Web: <http://www.clinisciences.com>



France

Company: CliniSciences S.A.S
Address: 74 Rue des Sulauses, 92000
Nanterre - France
Telephone: +33 9 77 40 09 09
Fax: +33 9 77 40 10 11
Email: info@clinisciences.com
Web: <http://www.clinisciences.com>



Switzerland

Company: CliniSciences Limited
Address: Markgasse 18 8302 Kloten -
Switzerland
Telephone: +41 (0)44 805 76 81
or +41 (0) 330 684 0982
Fax: +41 (0)1763 208 899
Email: switzerland@clinisciences.com
Web: <http://www.clinisciences.com>



UK

Company: CliniSciences Limited
Address: 11 Progress Business centre, White
Parkway, SL1 6DQ Slough - United Kingdom
Telephone: +44 (0)1753 866 611
or +44 (0) 330 684 0982
Fax: +44 (0)1763 208 899
Email: uk@clinisciences.com
Web: <http://www.clinisciences.com>



USA

Company: Biolrend Chemicals LLC
Address: c/o Carr Riggs Ingram,
600 Grand Boulevard, Suite 210 Miramar
Beach, FL 32650 - USA
Telephone: +1 860 650 7790
Fax: +1 860 650 4383
Email: info@biolrend-usa.com
Web: <http://www.biolrend-usa.com>



Iceland

Company: CliniSciences ApS
Address: Østertorvegade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: iceland@clinisciences.com
Web: <http://www.clinisciences.com>



Ireland

Company: CliniSciences Limited
Address: Ground Floor, 71 lower Baggot street
Dublin D02 P693 - Ireland
Telephone: +353 1 6971 146
Fax: +353 1 6971 147
Email: ireland@clinisciences.com
Web: <http://www.clinisciences.com>



Denmark

Company: CliniSciences ApS
Address: Østertorvegade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: denmark@clinisciences.com
Web: <http://www.clinisciences.com>



Germany

Company: Biolrend Chemistries GmbH
Address: Wilhelm-Meuser-Str. 4+43,
60927 Köln - Germany
Telephone: +49 221 9498 320
Fax: +49 221 9498 325
Email: info@biolrend.com
Web: <http://www.biolrend.com>



Italy

Company: CliniSciences S.r.l.
Address: Via Maremmana inferiore 378
Roma 00012 Guidonia Montecelio - Italy
Telephone: +39 06 94 80 56 71
Fax: +39 06 94 80 00 21
Email: italia@clinisciences.com
Web: <http://www.clinisciences.com>

